



ADVENTURER™

Semi-Micro, Analytical, and Precision Balances



Intuitive Balances Designed for Routine Weighing

Striking the ideal balance between inventive features and functional, uncomplicated weighing capabilities, the OHAUS Adventurer incorporates all of the applications necessary for routine weighing and measurement activities. With a color touchscreen, three level user management to fulfill GLP/GMP compliance capabilities, two USB ports, and much more, Adventurer is the most complete balance in its class.

Unique Features Include:

- Adventurer balances feature a color touchscreen, icon-based user interface, and an ergonomic design -making them easy to configure and use.
- Features such as specialized weighing modes, multiple connectivity options, and AutoCal™ provide versatility and flexibility for a variety of applications.
- Durable construction, large weighing surfaces, a space-saving draftshield design, and full housing in-use cover allow for use in lab, education and industrial environments.

ADVENTURER™ *Semi-Micro, Analytical, and Precision Balances*

Stability, Accuracy, and Fast Operation Ensure Optimal Weighing Results in Routine Weighing Tasks

Weighing Performance

- Delivers stable and reliable weighing results for routine weighing tasks

Stabilization Time

- Adventurer's fast stabilization time improves productivity in the laboratory

Calibration

- AutoCal™ — Selected models feature OHAUS' automatic internal calibration system that performs routine maintenance by calibrating the balance daily
- External Calibration — Traditional calibration in which the operator manually calibrates the balance with their choice of calibration weight value to ensure accuracy available on every model



Color Touchscreen Offers Easy and Fast Operation of Adventurer's Applications

- Operate and access Adventurer's nine application modes and abundant features that eliminate the need to do several manual calculations through the modern color touchscreen
- Operators can wear laboratory gloves while utilizing the touchscreen, eliminating the inconvenience and hazards associated with constantly putting on and removing gloves
- In addition to the touchscreen, Adventurer also has six mechanical keys that provide tactile feedback and allow the operator to perform repetitive operations such as tare, zero, calibration, and print



Application Modes



Weighing

Determine the weight of items in the selected unit of measure.



Parts Counting

Count samples of uniform weight.



Percent Weighing

Measure the weight of a sample displayed as a percentage of a pre-established Reference Weight.



Dynamic Weighing

Weigh an unstable load. Scale takes an average of weights over a period of time.



Density Determination

Determine density of solids or liquid. With the weigh below hook, it's possible to perform specific gravity tests for objects that cannot be easily placed on the weighing pan.



Check Weighing

Compare the weight of a sample against target limits.



Display Hold

Manually holds the last stable weight or highest weighing value on the display.



Totalization / Statistics

Measure cumulative weight of multiple items. Cumulative total may exceed balance capacity.



Formulation

For compounding and recipe making. The number of components can range from 2 to 50.



Batch Printing

Combine multiple samples into one printout rather than printing them one at a time.

ADVENTURER™ *Semi-Micro, Analytical, and Precision Balances*

Equipped with the Connectivity and Functional Features Required in Laboratories

Dual USB Ports

- A front USB host port is easily accessible and makes it simple to load data from the balance on to a flash drive without having to reach around to the back or move the balance
- A second USB device is located at the rear of the balance that can be used to connect the balance to a PC
- The connectivity options help meet traceability requirements in traditional installations

Label Printing Function

- Easy to link with Zebra printer and have one built-in label printing template

Balance Profiles

- The cloning feature allows you to save user and application settings to a USB flash drive which can be easily used to configure additional Adventurer balances

Below Minimum Sample Weight Indication

- When using the minimum weight feature, the display clearly indicates that your current sample weight is below your defined minimum limit. Simply increase your sample weight to assure that your results are up to your standard



Space-saving Draftshield Designed to Improve User Experience and Accessibility

- Draftshield doors are constructed of two glass panels, reducing the space required on the lab bench when the doors are open
- Wide door entry provides unobstructed access and allows larger weighing vessels to be easily placed on the pan, reducing the chance of accidental spillage
- Easy to keep clean in order to ensure a safe workspace by minimizing contamination



Power Saving Functions

- The Adventurer is designed with various power-saving features that help reduce the environmental impact. These include auto-off and brightness controls.



User Management Function ensures data security and data traceability

- 3 level user management function ensures data security and data traceability requirements
- One administrator, two supervisors and 10 users have preset accessibility in the software



Real Time Clock with GLP/GMP Data

- A real-time clock function keeps accurate time even during power loss
- GLP data capability has the ability to record Sample name, Project names and Balance IDs to help meet traceability and compliance requirements

ADVENTURER™

Semi-Micro, Analytical, and Precision Balances

Specifications

Model	AX125D	AX85	AX225D	AX124	AX224	AX324	AX223	AX423	AX523	AX623
External Calibration	-	-	-	AX124/E	AX224/E	-	AX223/E	AX423/E	AX523/E	AX623/E
Capacity (g)	82/120	82	102/220	120	220	320	220	420	520	620
Readability d, Fine Range (g)	0.00001			—			—			
Readability d, Full Range (g)	0.0001	0.00001	0.0001	0.0001			0.001			
Repeatability (sd.), ≤5% of Full Load (g)	0.00001			0.00008			0.0008			
Repeatability (sd.), 5% of Full Load to Fine Range Max (g)	0.00002			—						
Repeatability (sd.), Fine Range Max to Full Range (g)	0.0001	0.00002	0.0001				0.001			
Linearity Deviation, Typical (g)	±0.00006						±0.0006			
Linearity Deviation (g)	±0.0001			±0.0002			±0.002			
Stabilization Time (sec)	≤8			≤3			≤2			
Sensitivity Drift (PPM/°C)	2			3						
Min-Weight (Typical) (g) (USP, K=2, U=0.10%)	0.02			0.16			1.6			
Min-Weight (Optimal) (g) (USP, K=2, U=0.10%, SRP≤0.41d)*	0.0082			0.082			0.82			
Weighing Units	gram, milligram, mesgal, momme, Newton, ounce, pennyweight, Baht, carat, grain, pound,Tael (Hong Kong), Tael, (Singapore), Tael (Taiwan), tical, tola, troy ounce, custom (1)									
Weighing Applications	Weighing, Parts Counting, Percent Weighing, Check Weighing, Dynamic Weighing, Formulation, Density Determination, Totalization/Statistic, Display Hold, Batch Printing									
Pan Size	Ø 80 mm			Ø 90 mm			Ø 130 mm			
Calibration	All models feature external calibration. Models feature AutoCal™ internal calibration, except for AX.../E models									
Tare Range	To capacity by subtraction									
Power Requirements	AC Adapter Input: 100-240 VAC 0.3A 50-60 Hz AC Adapter Output: 12 VDC 0.84A									
Display Type	Full-Color Touchscreen WQVGA Graphic LCD									
Display Size	109 mm (diagonal)									
Base Housing (W×H×D)	354 × 230 × 340 mm									
Communication	RS232, USB Device, USB Host									
Temperature Range	10°C to 30°C									
Humidity Range	Maximum relative humidity 80 % for temperatures up to 30°C									
Storage Conditions	-10°C to 60°C at 10% to 90% relative humidity, non-condensing									
Shipping Dimensions	507 × 387 × 531 mm									
Net Weight	5.1 kg					5.8 kg				
Shipping Weight	7.8 kg					8.5 kg				


Model	AX622	AX1502	AX2202	AX4202	AX5202	AX6202	AX4201	AX8201	AX12001
External Calibration	AX622/E	AX1502/E	AX2202/E	AX4202/E	-	AX6202/E	AX4201/E	AX8201/E	AX12001/E
Capacity (g)	620	1520	2200	4200	5200	6200	4200	8200	12000
Readability d (g)	0.01						0.1		
Repeatability (sd.), ≤5% of Full Load (g)	0.008						0.08		
Repeatability (sd.), 5% of Full Load to Full Range (g)	0.01						0.1		
Linearity Deviation, Typical (g)	±0.006						±0.06		
Linearity Deviation (g)	±0.02						±0.2		
Stabilization Time (sec)	≤1.5								
Sensitivity Drift (PPM/°C)	3						5		
Min-Weight (Typical) (g) (USP, K=2, U=0.10%)	16						160		
Min-Weight (Optimal) (g) (USP, K=2, U=0.10%, SRP≤0.41d)*	8.2						82		
Weighing Units	gram, milligram, mesgal, momme, Newton, ounce, pennyweight, Baht, carat, grain, pound,Tael (Hong Kong), Tael, (Singapore), Tael (Taiwan), tical, tola, troy ounce, custom (1)								
Weighing Applications	Weighing, Parts Counting, Percent Weighing, Check Weighing, Dynamic Weighing, Formulation, Density Determination, Totalization/Statistic, Display Hold, Batch Printing								
Pan Size	175 x 195								
Calibration	All models feature external calibration. Models feature AutoCal™ internal calibration, except for AX.../E models								
Tare Range	To capacity by subtraction								
Power Requirements	AC Adapter Input: 100-240 VAC 0.3A 50-60 Hz AC Adapter Output: 12 VDC 0.84A								
Display Type	Full-Color Touchscreen WQVGA Graphic LCD								
Display Size	109 mm (diagonal)								
Base Housing (W×H×D)	354 × 230 × 100 mm								
Communication	RS232, USB Device, USB Host								
Temperature Range	10°C to 30°C								
Humidity Range	Maximum relative humidity 80 % for temperatures up to 30°C								
Storage Conditions	-10°C to 60°C at 10% to 90% relative humidity, non-condensing								
Shipping Dimensions	557 × 392 × 301 mm								
Net Weight	4.6 kg						3.8 kg		
Shipping Weight	6.5 kg						5.7 kg		

* The value for SRP is the standard deviation for n replicate weighings (n ≥ 10)


ADVENTURER™ *Semi-Micro, Analytical, and Precision Balances*

Outline Dimensions

0.1 and 1mg Models

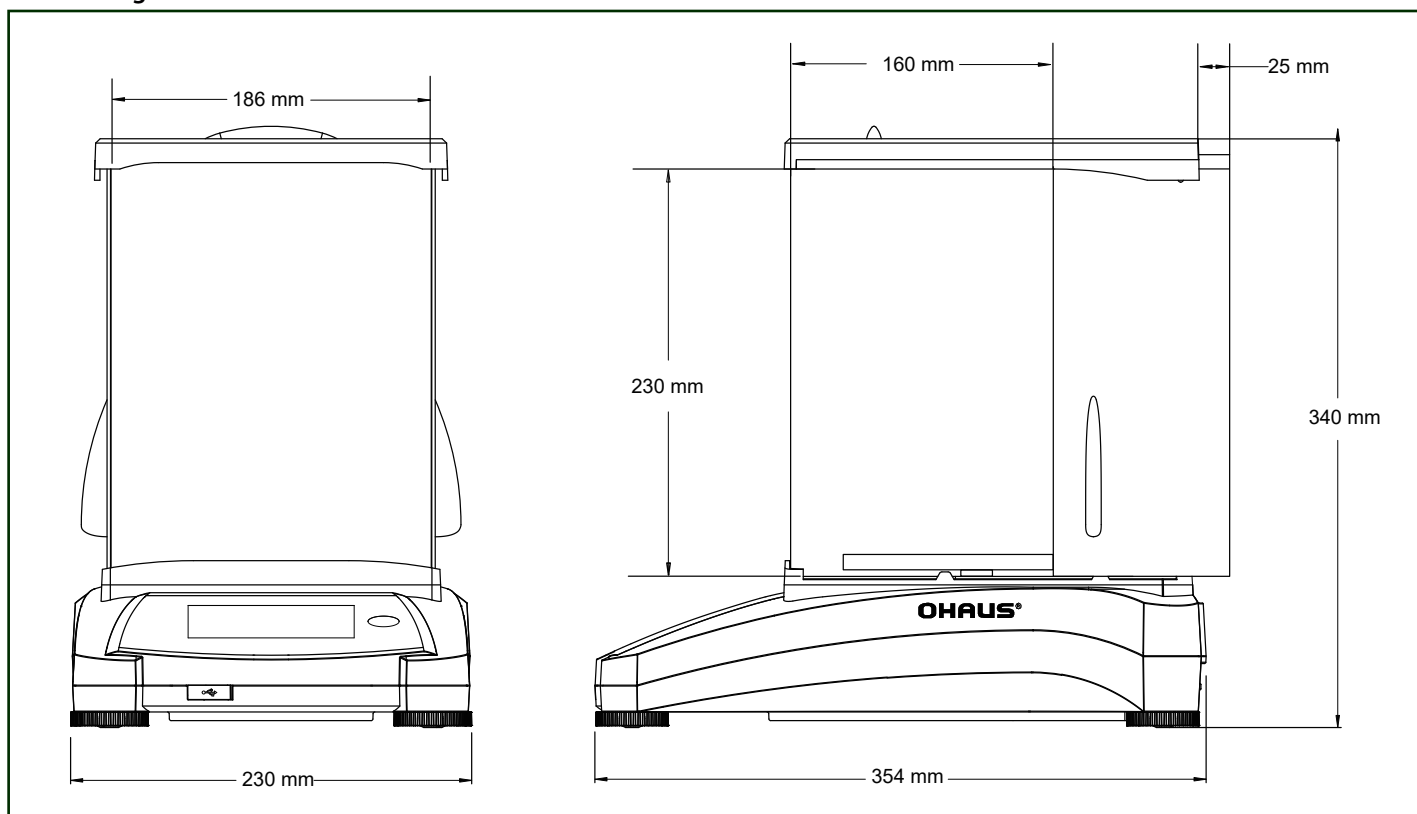

0.01 and 0.1g Models

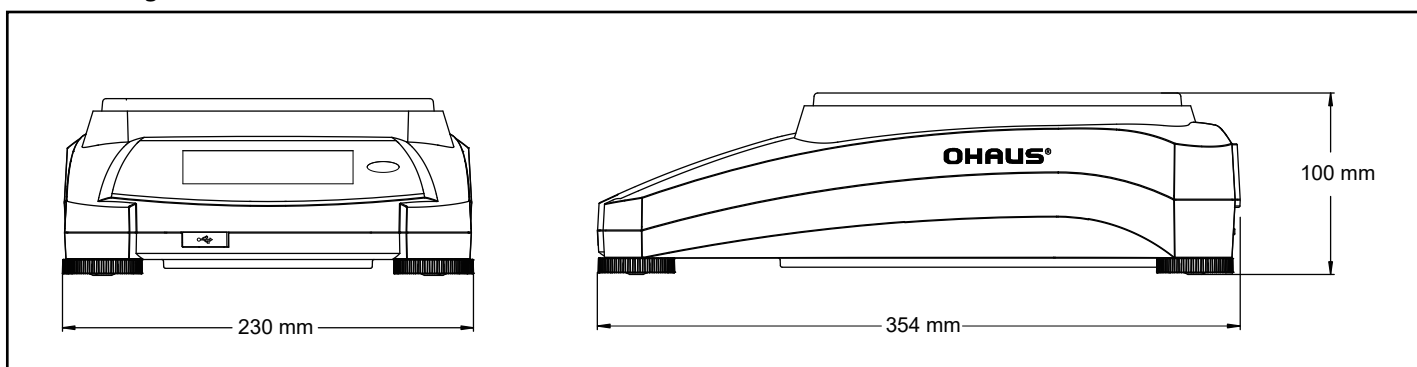

Additional Features

RS232 interface, integrated weigh below hook, full housing in-use cover, removable stainless steel pan, die-cast metal bottom housing, security bracket, illuminated up-front level indicator, four adjustable feet, software lockout menus, stability indicator, software overload/underload indicators, user selectable environmental settings, audible indicator, user selectable brightness settings, auto dim, auto-standby, auto-off, touchscreen calibration, auto tare, user selectable operating language (14), compatible interface command with MT-SICS and ST protocol

Compliance

Metrology: OIML R76, EN 45501 (Class I, nmax 320000; Class II, nmax 62000)

Product Safety: EN 61010-1, IEC 61010-1

Electromagnetic Compatibility: IEC 61326-1, EN 61326-1 (emissions Class B, immunity basic requirements)

Accessories

SF40A Impact Printer.....	30045641
Auxiliary Display, AD7-RS	30472064
Density Determination Kit.....	80253384
Cable, USB Interface (Type A to B).....	83021085
Security Device (Laptop Lock).....	80850043
RS232 Cable, PC 9 Pin	80500525
ION-100A AP Standalone Ionizer	30130303

OHAUS Asia/Pacific
Headquarters

e-mail: APmarketing@ohaus.com
Tel: +86 21 64855408

*With offices throughout
Europe, Asia, and
Latin America*

* ISO 9001:2008
Registered Quality
Management System

